

# VisiPro /VisiTrace DO Family

## Firmware Release History

### Firmware VisiTrace DO LDOUH006 and VisiPro DO ODOUH009

New Feature / Topics	Description	Advantage	Update
Bug fixed	Sensor data is lost from flash memory when powered by a noisy power supply	Sensor data is not lost from flash memory when powered by a noisy power supply	Recommended!

### Firmware VisiTrace DO LDOUH005 and VisiPro DO ODOUH008

New Feature / Topics	Description	Advantage	Update
Bug fixed	Reset of sensor, interrupt the 4-20 mA interface shortly and caused an error on the PCS (Process Control System)	Stable PCS communication	Recommended!

### Firmware VisiTrace DO LDOUH004 and VisiPro DO ODOUH007

New Feature / Topics	Description	Advantage	Update
Bug fixed	Flash memory loses the sensor data No S/N and Ref number are registered	No loss of sensor data	Recommended!

### Firmware LDOUH003 for VisiTrace DO only

New Feature / Topics	Description	Advantage	Update
Standardization of lifetime indication	Standardized lifetime indication	Reduce variation	Recommended!
Increase Product Calibration range from 20% to 50% of the reading	Now it is possible to perform product calibration with min./max. range of 50% of the reading		Recommended!

### Firmware ODOUH006 for VisiPro DO (Ex) only

New Feature / Topics	Description	Advantage	Update
Standardization of lifetime indication	Standardized lifetime indication	Reduce variation	Recommended!

## Firmware ODOUH005 for VisiPro DO (Ex) only

New Feature / Topics	Description	Advantage	Update
Fix Flash Memory Defective Error	Known problem when using the sensors in frequently power ups	No Flash Memory defective error	Recommended!
Bug fixes	Known bugs have been fixed		Recommended!

## Firmware LDOUH002 for VisiTrace DO only

New Feature / Topics	Description	Advantage	Update
Fix Flash Memory Defective Error	Known problem when using the sensors in frequently power ups	No Flash Memory defective error	Recommended!
Bug fixes	Known bugs have been fixed		Recommended!

© 2020 Hamilton Bonaduz AG. All rights reserved.

REF 695242/01 — 01/2020

# HAMILTON®

Web: [www.hamiltoncompany.com](http://www.hamiltoncompany.com)

USA: 800-648-5950

Europe: +41-58-610-10-10

### Hamilton Americas & Pacific Rim

4970 Energy Way  
Reno, Nevada 89502 USA  
Tel: +1-775-858-3000  
Fax: +1-775-856-7259  
[sales@hamiltoncompany.com](mailto:sales@hamiltoncompany.com)

### Hamilton Europe, Asia & Africa

Via Crusch 8  
CH-7402 Bonaduz, Switzerland  
Tel: +41-58-610-10-10  
Fax: +41-58-610-00-10  
[contact.pa.ch@hamilton.ch](mailto:contact.pa.ch@hamilton.ch)

To find a representative in your area, please visit [www.hamiltoncompany.com](http://www.hamiltoncompany.com).